

41-472100 ONE Broth Salmonella

*for the detection of Salmonella enterica with RiboFlow® Salmonella
Detection Kit within 24 hrs*

Components of ONE Broth Salmonella

1 x	ONE Broth Salmonella Base (500 g)	
4 x	Salmonella Additive ONE Broth	+2..+8 °C

Instructions for media preparation

- Dissolve 25 grams of ONE Broth Salmonella Base (Oxoid, CM1091) per litre of required medium in demineralized water.
- Autoclave medium at +121°C for 15 minutes and cool to below +50°C.
- Before use of the medium, dissolve the entire content of 1 vial of Salmonella Additive ONE Broth in 5 ml of sterile demineralized water.
- Per litre of medium, add 1 ml of dissolved additive.

Reconstituted additive is stable for 7 days at +2..+8°C

Reconstituted additive can be stored in aliquots at -20°C until the expiry date specified on the vial.