Instructions

Storage up
to 30 °C

Detection of coliforms in dairy products

Sample Preparation

Viable count in water or liquid foodstuff

1 mL of sample (diluted if necessary) is deposited on the CompactDry CF-L plate. Pre-treat the samples according to ISO 6887 or NMKL 91.

Viable count in solid food

Add buffer solution (e.g. 0.9 % saline solution, PBS, buffered peptone water) to the sample and homogenize using a Stomacher® (as per manufacturer instructions). Dispense 1 mL of sample (diluted if necessary) in the centre on the CompactDry CF-L plate.

Viable count in swab test sample

Use the swab to wipe the surface, put the swab back into the tube and screw tightly. Shake the tube to mix. Dispense 1 mL of sample (diluted if necessary) in the centre on the CompactDry CF-L plate.

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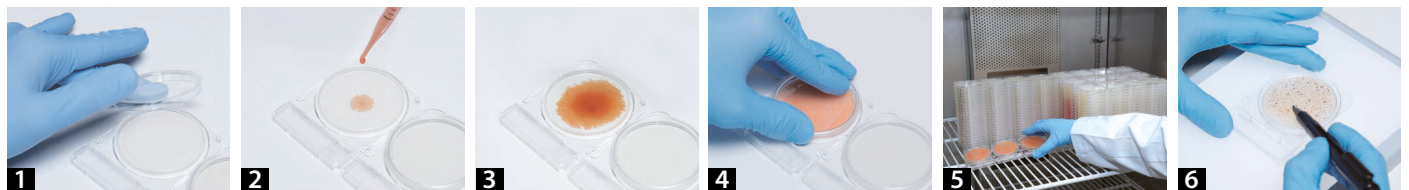
CompactDry Protocol

1 Open the cap **2** Dispense 1 mL of sample in the middle of the CompactDry plate OR add filter as instructed **3** Sample diffuses passively and evenly across the dehydrated media sheet transforming it into a gel within seconds **4** Replace the cap, label and turn the plate over **5** Incubate for the required time and temperature. **6** Following incubation, count the number of coloured colonies.

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Incubation temperature: 37 ±1 °C
Incubation time: 24 ±2 hours

Interpretation

Colonies formed are blue/blue green in colour.

General information

Growth of none coliforms bacteria is generally inhibited.

The growth area is 20 cm². The back of the plate has a 1 cm² grid to ease colony counting. If there are a large number of colonies, the total viable count can be obtained by multiplying the average number of colonies in several grids by 20.

High concentrations (>250 CF-LU) will cause the entire growth area to become white/light blue. In this case dilute the sample.

After use please follow local disposal regulations.

Storage and shelf life

Store at room temperature (+1 to +30 °C). Shelf life 24 months after manufacturing.

PACKAGES	PRODUCT NUMBER
40 plates	54065-CFL-0040
240 plates	54065-CFL-0240

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